

Add a Filter Wheel and a Shutter

Launch HCIImage, go to File and select Current Profile. In the Device Control tab, select Filter Devices, click Add and follow the instructions below.

The figure consists of four screenshots of the 'Add FILTER Device' dialog box, each showing a different tab and configuration step. Red dotted lines and numbered callouts (1-4) point to specific elements in each screenshot.

1 Device Control
Select the controller, the type of filter wheel and enable Add Shutter

2 Enable COM Port
Select the COM Port for the device

3 Filter Setup
Select the filter position, enter the name and click Modify. Repeat for each filter position

4 Shutter Setup
Set the status to Active and click OK

The dialog box has four tabs: Controller, COM-Port, Filter Setup, and Shutter Setup. The first screenshot shows the Controller tab with a list of controllers and filter wheel options. The second screenshot shows the COM-Port tab with a list of COM ports and communication settings. The third screenshot shows the Filter Setup tab with a list of filter positions and a 'Customize Positions' section. The fourth screenshot shows the Shutter Setup tab with a list of shutters and shutter settings.